

Macro-imager footpedal installation instructions v2.0

PLEASE NOTE: You must first upgrade the Macro-imager software to Version 2v22 in order to use the new footpedal. Please see Macro-imager software upgrade instructions v2.02 for more info.

1. Software downloads:

- Visit the Macro-imager website to download the software:
- <https://www.macro-imager.com/downloads.html>
- Download "Pedable" - software to program and control the footpedal
- Once downloaded, you will need to "unzip" it before running - pedablesetup.exe and piehid.exe

2. Installation - Footpedal:

Pedable:

1. Plug the footpedal into an available USB port on the PC.
2. Run pedablesetup.exe
3. Agree to the license then click OK (even if it does not detect the pedal).
4. Go to OPTIONS - select the **VEC Infinity USB Pedals**
5. You can select / deselect run on start-up. You can manually run Pedable as and when the footpedal is required if you prefer.

Pedable will automatically download an update to v2.01. If no Internet is available or if security software prevents this please:

RIGHT-CLICK on piehid.exe

Select "Run as administrator"

- Double-click the "Center" pedal (**CAPTURE**)
- Select SEND KEY COMBINATION - then Next
- Press CTRL then Q
- FINISH

Repeat as follows:



"Left" = CTRL and E (Continuous Zoom OUT)

"Right" = CTRL and W (Continuous Zoom IN)

OR - For CALIBRATION MODULE:

"Left" = CTRL and D (Step Zoom OUT)

"Right" = CTRL and S (Step Zoom IN)

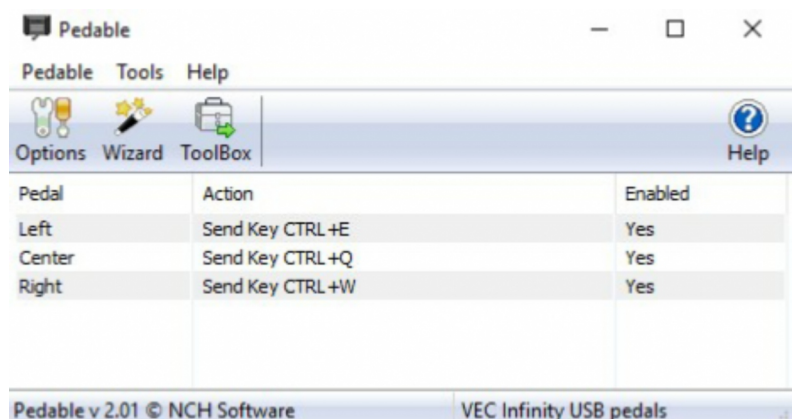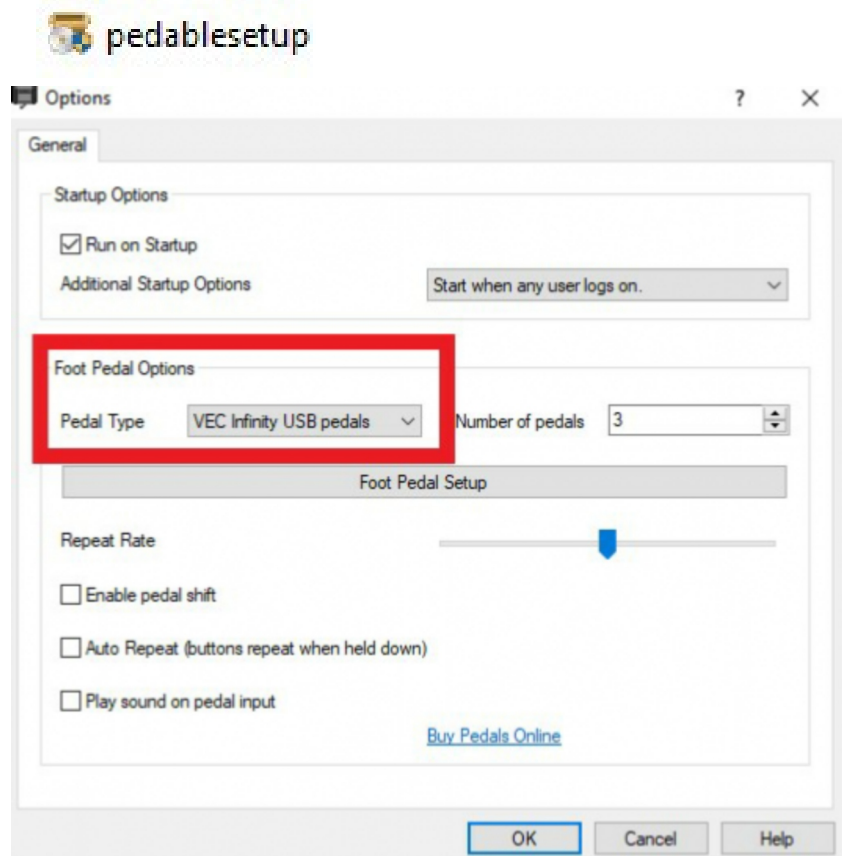

- Select the Pedable menu and Exit (close to tray) to finish.