

Macro-imager Upgrade Instructions v2.02

Before you begin the upgrade, please save a copy of the file SessionOptions.xml located in C:\Program Files (x86)\Micropix Ltd\Macro-imager as a back-up (e.g. on the desktop) - you will need it later.

1. Software downloads:

- Visit the Macro-imager website to download the software:
- <https://www.macro-imager.com/downloads.html>
- Download the latest Macro-imager software - this is a 32-bit application regardless of your operating system.

Once downloaded, you will need to “unzip” the software and save it.

2. Installation - Macro-imager:

- Go to the CONTROL PANEL and uninstall Macro-imager.

1. Open the MacroImagerInstaller_2v02 folder.
2. Double-click “MacroImagerInstaller” (**NOT SET-UP**).
3. Click “Next” 3 times and let the installer run.
4. When you get the **Installation Complete** message, click Close to finish.
5. Copy / cut your saved SessionOptions.xml file and put it back into C:\Program Files (x86)\Micropix Ltd\Macro-imager - please replace the file that is already there.

